


# D-Luciferin, Sodium Salt

## Product Sheet and Protocol

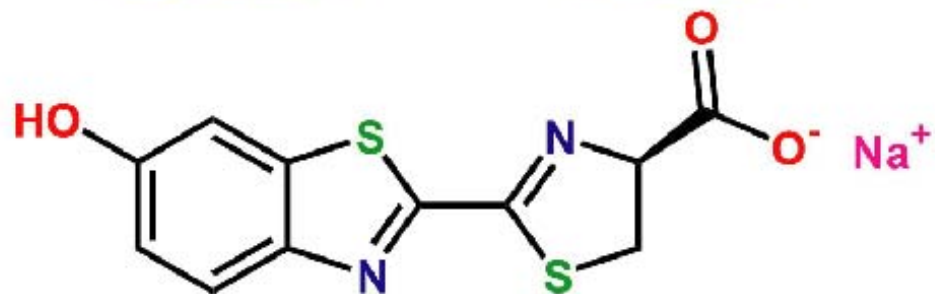

# D-Luciferin, Sodium Salt

## Instruction Manual

**D-Luciferin Sodium salt (Na<sup>+</sup>)** is designed for use in *in vitro* and *in vivo* bioluminescent assays. The quality and purity of the D-Luciferin is essential to obtain good and reproducible results. OZ Biosciences offers high quality, **Endotoxin-Free** D-Luciferin Na<sup>+</sup> salt.

### List of D-Luciferin products:

Catalog Number	Description	Weight
LK10000	D-Luciferin Potassium salt	1 g
LN10000	D-Luciferin Sodium salt	1 g

Use the table above to determine the appropriate catalogue number for your needs. You can order these products by contacting us. These products are also available in bigger quantities (2g, 5g, 10g etc.). Please contact us for a quotation at [order@ozbiosciences.com](mailto:order@ozbiosciences.com). For all other supplementary information, do not hesitate to contact our dedicated technical support ([tech@ozbiosciences.com](mailto:tech@ozbiosciences.com)).

#### OZ BIOSCIENCES

Parc Scientifique de Luminy  
Zone Luminy Entreprise  
163, avenue de Luminy – Case 922  
13288 Marseille Cedex 9 – France  
Tel: +33 (0)4.86.94.85.16  
Fax: +33 (0)4.86.94.85.15  
E-mail: [contact@ozbiosciences.com](mailto:contact@ozbiosciences.com)  
Web Site: [www.ozbiosciences.com](http://www.ozbiosciences.com)



**OZ BIOSCIENCES**  
The art of delivery systems

## 1. Description

### 1.1. Introduction and characteristics

Congratulations on your purchase of the **D-Luciferin, Sodium Salt** product!


**D-Luciferin, Sodium Salt** is a synthetic Firefly Luciferin. This highly pure and endotoxin-free product is perfect for *in vitro* and *in vivo* bioluminescent assays. The highest quality of substrate eliminates possible interference in assays due to the presence of endotoxins. Proper packaging in amber vials and under neutral gas assures product integrity and stability. This water soluble substrate of Luciferase is ready and easy to use.

#### **Main Features are:**

- High purity > 99.7%
- Good solubility and great sensitivity
- Reliable *in vivo* reporter for bioluminescent assays
- Endotoxin free (ideal for *in vivo* application)
- Suitable for *in vitro* experiments
- Easy to use
- Quick and easy distribution throughout the animal

#### **D-Luciferin, Sodium Salt:**

- *Chemical Name:* 4,5-Dihydro-2-(6-hydroxy-2-benzothiazolyl)-4-thiazolecarboxylic acid sodium salt
- *Molecular Formula:* C<sub>11</sub>H<sub>7</sub>N<sub>2</sub>O<sub>3</sub>S<sub>2</sub>·Na·H<sub>2</sub>O
- *Molecular weight:* 320.32 g/mol
- *CAS Number:* 103404-75-7
- *Molecular Structure:*

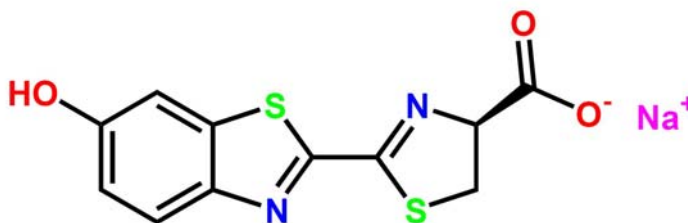

The molecular structures of Firefly and Beetle Luciferin are identical.

- *Molecular biology grade and premium pure:* 99.7% pure. Quality verified by nine independent criteria including HPLC and FTIR.

### 1.2. Kit content

D-Luciferin, sodium salt is provided in a 1.0g/vial in an amber bottle under nitrogen. D-Luciferin is sensitive to light, oxygen and moisture. This product is also available in larger quantities (2g, 5g, 10g etc.). Please contact us for a quotation at [order@ozbiosciences.com](mailto:order@ozbiosciences.com)

### 1.3. Stability and Storage

D-Luciferin is sensitive, to light, oxygen and moisture as powder and in solution.

#### **Storage (powder)**

Upon receipt and for long-term use, store the powder in a tightly closed and desiccated container at -20°C. D-Luciferin is stable for at least one year at -20°C under desiccated and neutral gas conditions. If the bottle containing the D-Luciferin powder has to be opened and closed repetitively, it is recommended to fill the vial with an inert gas (Argon or Nitrogen) before closing.

#### **Storage (solution)**

D-Luciferin stock solution can be prepared in water or D-PBS and then frozen at -20°C or preferably at -80°C. Stability of the D-Luciferin solution is rather controversial and accordingly we recommend for the most sensitive studies (low concentrations of enzyme, sub-optimal temperatures or salt concentrations) to use freshly made solutions. Nonetheless, frozen D-Luciferin solutions (-80°C) have been suitable for use after at least 6 months. D-Luciferin is quite unstable at low pH (< 6.5) and high pH (> 7.5). To avoid repeated cycles of freezing and thawing, we recommend aliquoting D-Luciferin solution in small tubes and storing them at -80°C.

#### **Shipping condition**

D-Luciferin is shipped at room temperature.

## 2. Applications

- Bioluminescent assays in living cells, tissues and animal models
- Luciferase reporter gene assays
- Whole animal imaging (*in vivo* reporter assay)
- Appropriate read-out for transfection/transduction with luciferase reporter gene and luciferase-fusion constructs
- ATP assays (Luciferase catalyzes conversion of ATP into AMP) and immunoassays
- Pyrosequencing
- Luciferase fragment complementation for sequential gene analysis experiments

## 3. Protocol

The instructions given below represent successfully applied protocols. They can be used as guidelines to quickly achieve very high bioluminescence signal. Optimal conditions do vary according to animals, cell cultures, route of administration and assay sensitivity. D-Luciferin final quantity might have to be adjusted to achieve best results.

### 3.1. Preparation of D-Luciferin for *in vivo* assay

Biodistribution of the D-Luciferin is rapid and easy throughout the animal but kinetics may be tissue dependent. D-Luciferin can penetrate cell membranes and is able to pass blood-brain barrier, blood-placenta barrier and blood-testis barrier.

#### **Materials**

- D-Luciferin, sodium salt (#LN10000)
- D-PBS (without Mg<sup>2+</sup> and Ca<sup>2+</sup>)
- Syringe filter 0.2 µM

#### **Procedure**

The best is to reconstitute the quantity of needed D-Luciferin for each experiment. However, stocks of frozen solution of Luciferin can also be used (*see storage and stability section above*).

1. Dissolve the D-Luciferin in D-PBS to obtain a final concentration of 15 or 30mg/mL.
2. Filter sterilize through a 0.2µM filter (optional, not required).
3. Inject intraperitoneally or intravenously 5 to 15 minutes before imaging each animal with 10µL/g of body weight or 150 to 300 µg/g of body weight. To have a sufficient or even excess amount of substrate we recommend using 3mg of D-Luciferin per mice (100µL of 30mg/mL).

**Note 1:** we recommend performing *kinetic study* to determinate D-Luciferin kinetic curve and peak signal for your animal model:

- i. Inject D-Luciferin as previously described. Awake or sedated animals can be used; however, with sedated animals the kinetics (peak luciferase expression time) may slightly be extended.
- ii. After 3 min. if awoken animals were used in step 1, proceed to sedate animals.
- iii. Place sedated animals in imaging chamber and capture the first image about 5 min. after the Luciferin injection.
- iv. Continue to take images every 5-10 min. up to about 45-60 min. to produce a kinetic curve for Luciferin expression in your animal model.
- v. Thereafter you can choose the best time point to image at. We image most of our models at 10-20 min. after D-Luciferin injection.

**Note 2:** for Intraperitoneal (I.P.) injection of D-Luciferin.

The animal needs to be in dorsal position (abdomen side up), manually controlled, with cranial end of animal pointed down. Injection has to be made in the animal's lower left abdominal quadrant. Needle (25 gauge) should be bevel-side up and slightly angled when entering the abdominal cavity. Penetrate just through abdominal wall (about 4-5 mm) and inject (1cc syringe).

### 3.2. Preparation of D-Luciferin for *in vitro* assay

#### Material

- D-Luciferin, sodium salt (#LN10000)
- Molecular biology grade water
- Complete media

#### Procedure

1. Prepare a D-Luciferin stock solution of 15 or 30mg/mL (100X or 200X) in molecular biology grade water
2. Mix by inverting the tube or bottle repeatedly until D-Luciferin is completely dissolved
3. At this step, solution can be immediately used or aliquoted and frozen at -20°C or -80°C
4. Add stock solution of D-Luciferin to pre-warmed tissue culture medium, the final D-Luciferin concentration should be 150µg/mL.
5. Remove old cell culture medium from the cultured cells
6. Add the D-Luciferin solution (150µg/mL) to cells immediately before imaging

*Note: Cells can be incubated at 37°C for a short period of time before imaging to increase the signal*

## 4. Related Products

Description	Reference
<b>Magnetofection Technology</b>	
Mega Magnetic Plate	MF14000
Super Magnetic Plate	MF10000
PolyMag 1mL ( <i>for all nucleic acids</i> )	PN31000
PolyMag Neo 1mL ( <i>for all nucleic acids</i> )	PG61000
CombiMag 1mL ( <i>to boost transfection reagent</i> )	CM21000
SilenceMag 1mL ( <i>for siRNA application</i> )	SM11000
NeuroMag 1mL ( <i>for transfection of neurons</i> )	NM51000
ViroMag 1mL ( <i>for all viral applications</i> )	VM41000
ViroMag R/L 1mL ( <i>for retrovirus and lentivirus</i> )	RL41000
AdenoMag 1mL ( <i>for adenovirus</i> )	AM71000
<b>Protein Delivery Systems</b>	
Ab-DeliverIN 1 mL	AI21000
Pro-DeliverIN 1 mL	PI11000
<b>Lipofection (lipid-based reagents)</b>	
Lullaby siRNA transfection reagent	LL71000
DreamFect Gold Transfection reagent 1mL	DG81000
DreamFect Transfection reagent 1mL	DF41000
VeroFect Transfection Reagent 1mL	VF61000
FlyFectin Transfection Reagent 1mL	FF51000
<b>CaPO Transfection Kit</b>	CP90000
<b>Plasmids pVectOZ</b>	
pVectOZ-LacZ 25µg	PL00030
pVectOZ-Luc 25µg	PL00040
pVectOZ-GFP 25µg	PL00020
<b>Gene &amp; Protein Tools</b>	
Bradford – Protein Assay Kit	BA00100
β-Galactosidase (ONPG) assay kits	GO10001
β-Galactosidase (CPRG) assay kits	GC10002
X-Gal Staining Kit	GX10003

Our dedicated and specialized technical support group will be pleased to answer any of your request and to assist you in your experiments. Do not hesitate to contact us for all complementary information and remember to visit our website in order to stay inform on our last breakthrough technologies and updated on our complete product list. <http://www.ozbiosciences.com>.

## Purchaser Notification

### Limited License

The purchase of the **D-Luciferin, sodium salt** product grants the purchaser a non-transferable, non-exclusive license to use the kit and/or its separate and included components (as listed in section 1, Kit Contents). This reagent is intended **for internal research only** by the buyer. Such use is limited to the use described in the product manual. In addition, research only use means that this kit and all of its contents are excluded, without limitation, from resale, repackaging, or use for the making or selling of any commercial product or service without the written approval of OZ Biosciences.

Separate licenses are available from OZ Biosciences for the express purpose of non-research use or applications of the **D-Luciferin, sodium salt** product. To inquire about such licenses, or to obtain authorization to transfer or use the enclosed material, contact the Director of Business Development at OZ Biosciences.

Purchasers may end this License at any time by returning all **D-Luciferin, sodium salt** product and documentation to OZ Biosciences, or by destroying all **D-Luciferin, sodium salt** product. Purchasers are advised to contact OZ Biosciences with the notification that a **D-Luciferin, sodium salt** product is being returned in order to obtain a refund and/or to expressly terminate a research only license granted through the purchase of the kit(s).

This document covers entirely the terms of the **D-Luciferin, sodium salt** product only license, and does not grant any other express or implied license. The laws of the French Government shall govern the interpretation and enforcement of the terms of this License.

### Product Use Limitations

The **D-Luciferin, sodium salt** product and all of its components are developed, designed, intended, and sold for research use only. They are not to be used for human diagnostic or included/used in any drug intended for human use. All care and attention should be exercised in the handling of the kit components by following proper research laboratory practices.

For more information, or for any comments on the terms and conditions of this License, please contact:

Director of Business Development  
OZ Biosciences  
Parc Scientifique de Luminy  
Zone Luminy Entreprise  
163, avenue de Luminy – Case 922  
13288 Marseille Cedex 9 - FRANCE  
Tel: +33 (0)4.86.94.85.16  
Fax: +33 (0)4.86.94.85.15  
E-mail: [business@ozbiosciences.com](mailto:business@ozbiosciences.com)